

**Instruction Sheet
for the PASCO
Model SE-9415**

4mm BANANA PATCHCORD KIT

Introduction


This patch cord kit has been supplied in pieces so that one may custom design a patch cord to any length. Use the setup procedure to assemble parts.

Additional Equipment Required:


- Screw driver (narrow, standard blade)
- Wire cutter
- Wire stripper
- Soldering equipment
- Emery cloth, sandpaper, or small file

Setup Procedure


- ① Cut wire to desired length. Strip 1/4" of wire jacket from each end of wire.

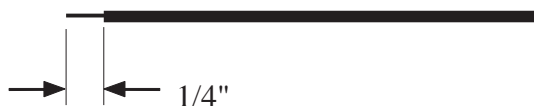


- ② Use the narrow screwdriver to pry open the end cover. Remove banana plug from plug housing by opening end cover and pushing banana plug through hole in housing.

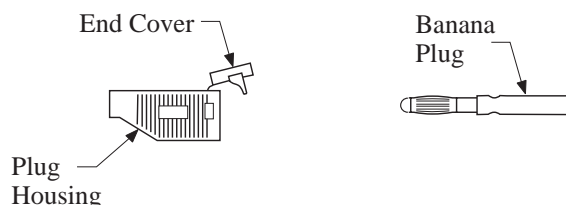

- ③ Use emery cloth, fine sandpaper or a file to roughen the surface of the banana plug near one of the side holes so the solder will stick. Place exposed wire into large thru-hole in banana plug and solder.

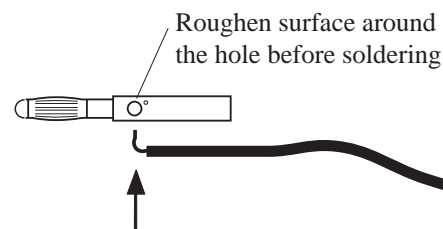

- ④ While holding banana plug and wire together, insert through hole in plug housing until plug seats in hole.

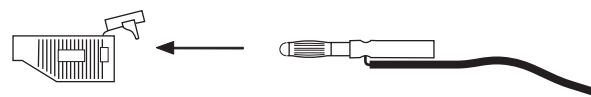

- ⑤ Align wire with hole in end cover and snap end cover closed.



Limited Warranty

PASCO scientific warrants this product to be free from defects in materials and workmanship for a period of one year from the date of shipment to the customer. PASCO will repair or replace, at its option, any part of the product which is deemed to be defective in material or workmanship. This warranty does not cover damage to the product caused by abuse or improper use. Determination of whether a product failure is the result of a manufacturing defect or improper use by the customer shall be made solely by PASCO scientific. Responsibility for the return of equipment for warranty repair belongs to the customer. Equipment must be properly packed to prevent damage and shipped postage or freight prepaid. (Damage caused by improper packing of the equipment for return shipment will not be covered by the warranty.) Shipping costs for returning the equipment, after repair, will be paid by PASCO scientific.

Limited Warranty

PASCO scientific warrants this product to be free from defects in materials and workmanship for a period of one year from the date of shipment to the customer. PASCO will repair or replace, at its option, any part of the product which is deemed to be defective in material or workmanship. This warranty does not cover damage to the product caused by abuse or improper use. Determination of whether a product failure is the result of a manufacturing defect or improper use by the customer shall be made solely by PASCO scientific. Responsibility for the return of equipment for warranty repair belongs to the customer. Equipment must be properly packed to prevent damage and shipped postage or freight prepaid. (Damage caused by improper packing of the equipment for return shipment will not be covered by the warranty.) Shipping costs for returning the equipment, after repair, will be paid by PASCO scientific.