




Basic Dynamics Systems

ME-5701, ME-5702, ME-5703, ME-5704, ME-5705, ME-5706

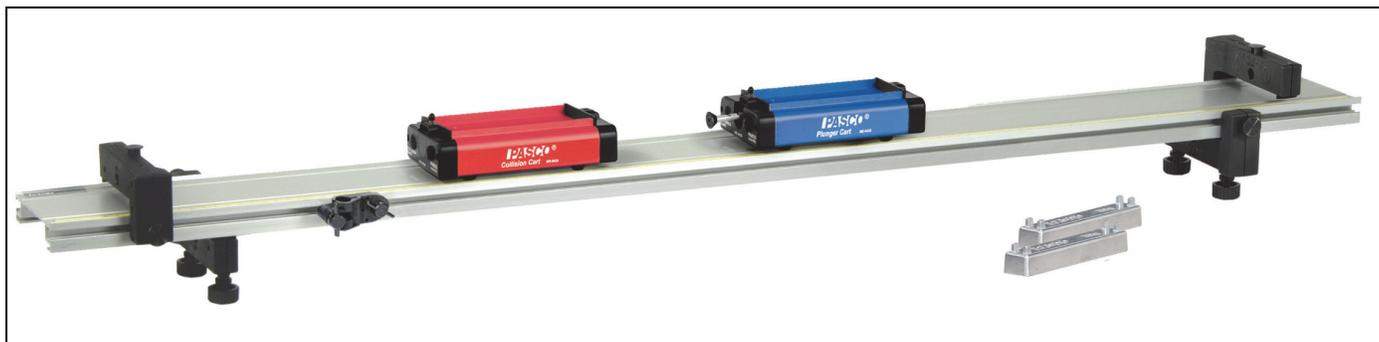

Basic Systems	Plastic Track	1.2 m Metal Track	2.2 m Metal Track
Plastic Carts	ME-5701	ME-5702	ME-5703
Metal Carts	ME-5704	ME-5705	ME-5706


Introduction

The Basic Dynamics Systems vary by the type of track and carts they include. They have in common that a variety of experiments and demonstrations can be done. You can also combine them with many other PASCO products—both traditional and computer—based—for an even greater range of uses.

This manual contains descriptions of the included equipment, complete instructions for five experiments, and several additional experiment suggestions.


Included Equipment

ME-5701	Items
ME-6757A	2 Masses (250 g)
ME-6950	PAScar (set of 2)
ME-6960	PAstrack
ME-8971	End Stops (set of 2)
ME-9836	Pivot Clamp

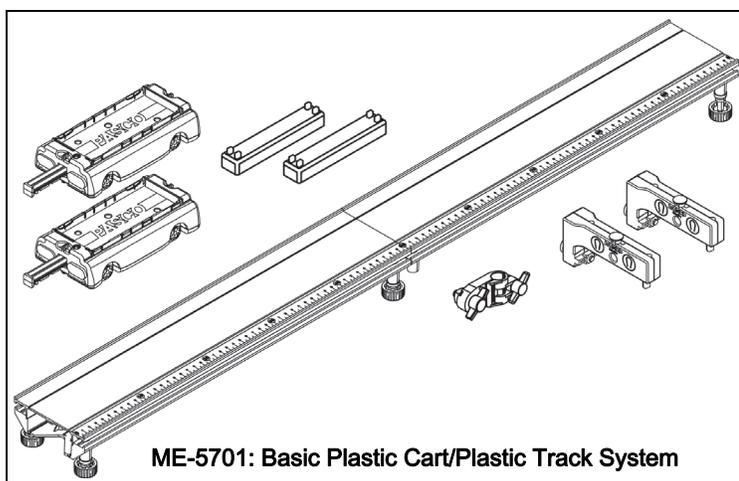


ME-5702 Items

ME-6757A	2 Masses (250 g)
ME-6950	PAScar (set of 2)
ME-8971	End Stops (set of 2)
ME-8972	Adjustable Feet (2)
ME-9493	Metal Track 1.2 m
ME-9836	Pivot Clamp

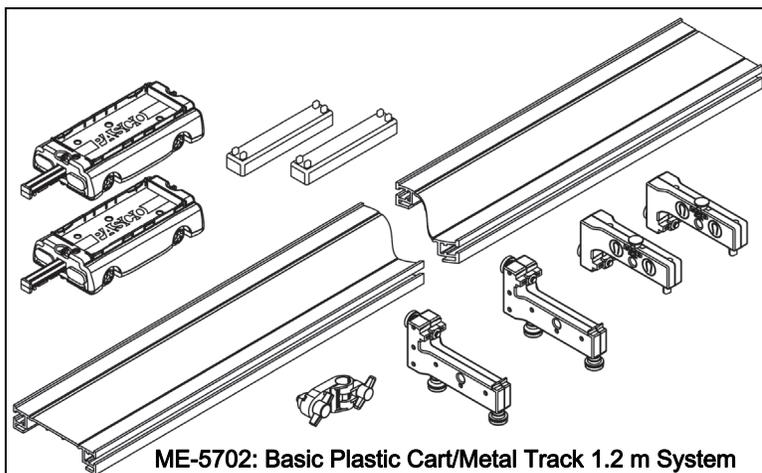
**ME-5702: Basic Plastic Cart/Metal Track 1.2 m System****ME-5703 Items**

ME-6757A	2 Masses (250 g)
ME-6950	PAScar (set of 2)
ME-8971	End Stops (set of 2)
ME-8972	Adjustable Feet (2)
ME-9779	Metal Track 2.2 m
ME-9836	Pivot Clamp

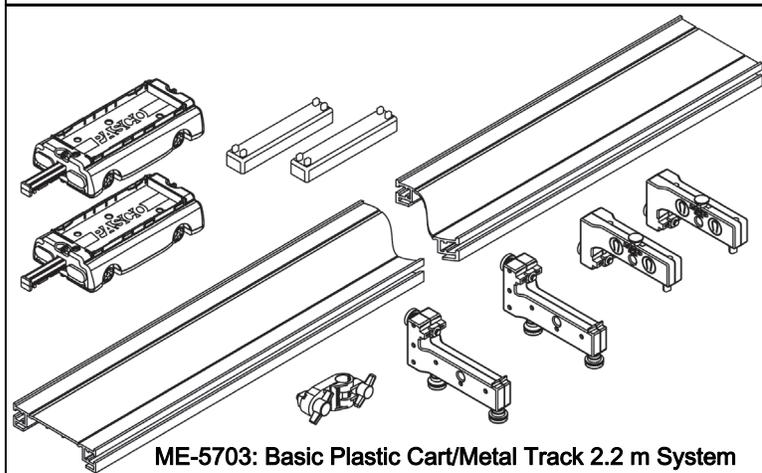**ME-5703: Basic Plastic Cart/Metal Track 2.2 m System****ME-5704 Items**

ME-6757A*	4 Masses (250 g)
ME-6960	PAStreck
ME-8971	End Stops (set of 2)
ME-9430	Plunger Cart
ME-9454	Collision Cart
ME-9836	Pivot Clamp


*Quantity 2

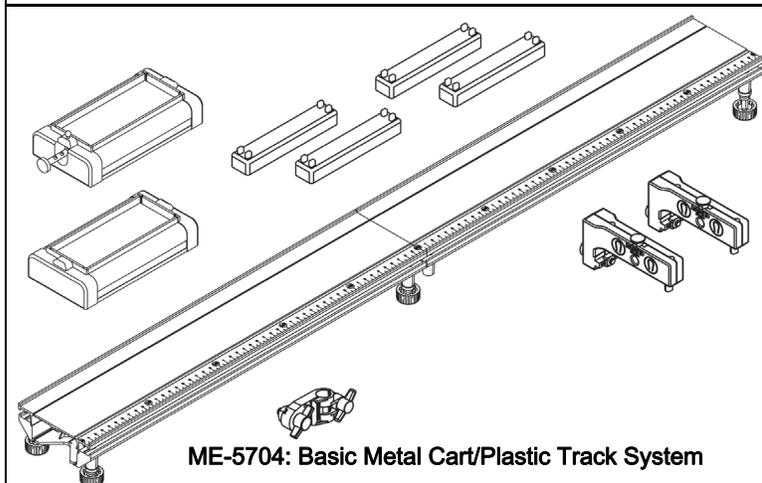**ME-5704: Basic Metal Cart/Plastic Track System**

ME-5705 Items

- ME-6757A* 4 Masses (250 g)
- ME-8971 End Stops (set of 2)
- ME-8972 Adjustable Feet (2)
- ME-9430 Plunger Cart
- ME-9454 Collision Cart
- ME-9493 Metal Track 1.2 m
- ME-9836 Pivot Clamp


*Quantity 2

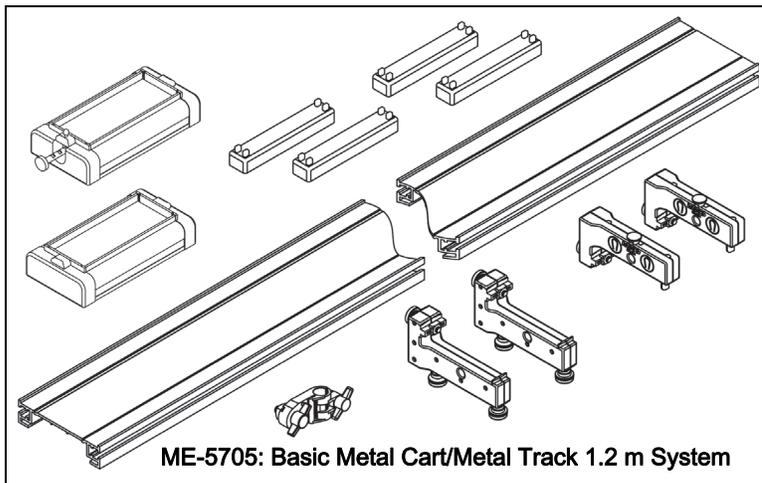

ME-5706 Items

- ME-6757A* 4 Masses (250 g)
- ME-8971 End Stops (set of 2)
- ME-8972 Adjustable Feet (2)
- ME-9430 Plunger Cart
- ME-9454 Collision Cart
- ME-9779 Metal Track 2.2 m
- ME-9836 Pivot Clamp

*Quantity 2

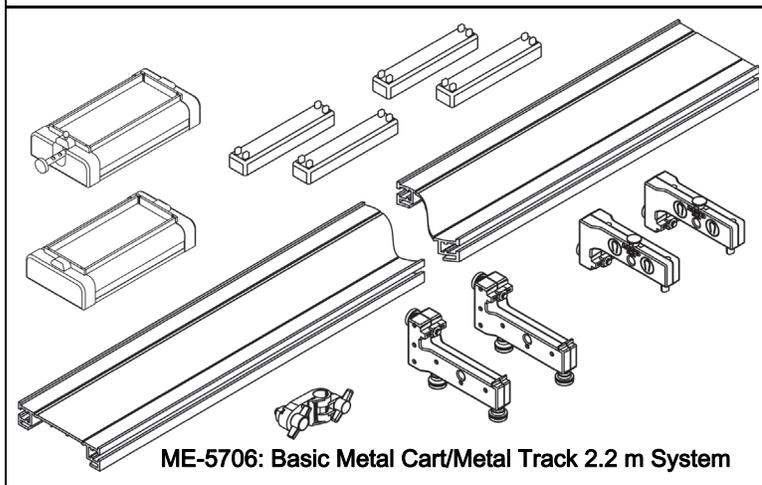


About the Equipment

Metal Track (1.2 m or 2.2 m) in ME-5702, ME-5703, ME-5705 and ME-5706

Metal Track The aluminum track has two grooves to guide the wheels of carts, a metric scale for measuring cart positions, and T-slots on both sides for attaching end stops, leveling feet and other accessories.

Tip: The track is designed to support the weight of carts and related apparatus. Excess weight will warp it. When you store the track, ensure that no heavy object will be placed on top of it.

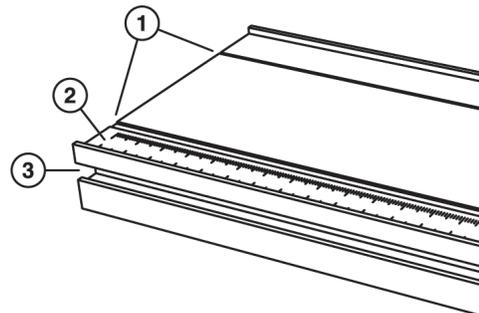



Track Features

1: Grooves 2: Metric scale 3: T-slot

Adjustable Feet

Attach the feet to the track as illustrated. Turn the feet screws to level the track then tighten the lock nuts to secure them.

For maximum stability, position the feet about 1/4 of the track length from each end

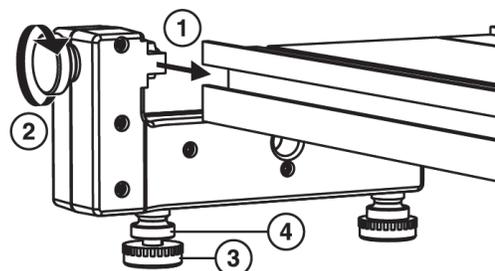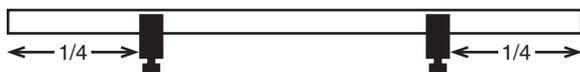


Adjustable Feet

- 1: Slide tab into T-slot.
- 2: Tighten thumbscrew to secure feet set.
- 3: Turn feet screws to level track.
- 4: Tighten lock nuts to secure feet screws.

Adjustable End Stops

Attach the end stops to the track as illustrated. The end stops contain magnets to repel the magnetic bumper of a cart. A stud is provided as an attachment point for a spring.

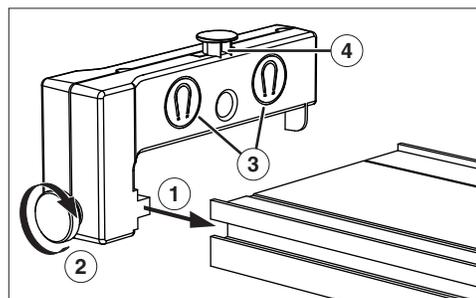

Adjustable End Stop

- 1: Slide tab into T-slot on side of track.
- 2: Tighten thumbscrew to secure end stop.
- 3: Magnets repel cart's magnetic bumper
- 4: Spring attachment stud

Plastic Track in ME-5701 and ME-5704

PAStack The PAStack consists of two sections that can be joined together with the included connector and split apart for storage. Together the two sections form a 1 m track. An extra connector is included, allowing any number of tracks to be joined together.

To assemble the track:

1. Slide a connector piece into the “50 cm” end of one track section. Push it hard until you hear it snap.
2. Slide the other track section onto the connector piece. Push it until it snaps.
3. Turn the track upside down and tighten the screws on the connector.


To level the track:

1. Set the track upright on a level surface.
2. Place a cart on the track to test it for level.
3. Adjust the feet to make the track straight and level.
4. Tighten the locknut of each foot.

To disassemble the track:

1. Loosen the connector screws.
2. Press and hold the release tabs on the bottom of the track.
3. Pull the track sections apart.

For more information, see the instruction sheet packaged with the track.

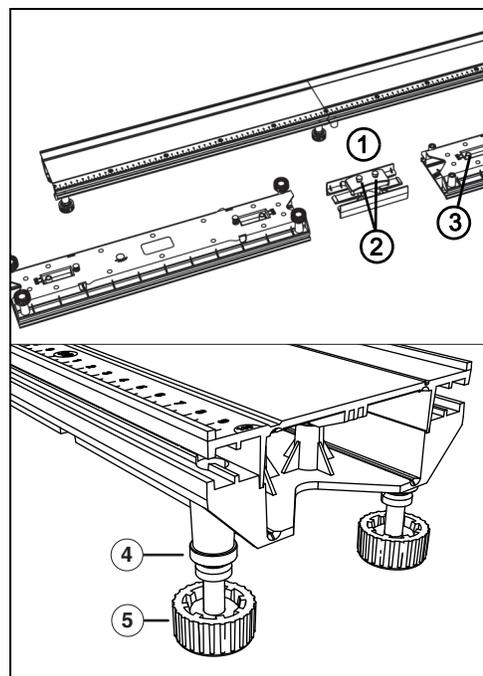

PAStack Parts

- 1: Connector
- 2: Connector screws
- 3: Release tabs
- 4: Lock nut
- 5: Foot

Plastic Carts in ME-5701, ME-5702, and ME-5703

PAScars These carts run along the track on low-friction wheels.


Each cart includes a spring-loaded plunger that launches the cart off of an end stop or another cart. Push the plunger in to the first, second, or third “click” (for a slow, medium, or fast launch). To make the plunger pop out, press the trigger button. When the plunger is not in use, push it all the way in to the fourth position for storage.

For elastic and inelastic collisions, both carts have hook-and-loop bumpers on one end and magnetic bumpers on the other end. When the hook-and-loop bumpers collide, the carts stick to each other. When the magnetic bumpers collide, the carts repel each other without touching. The magnetic bumpers also repel the magnets in the end stops.

The top of each cart forms a tray designed to hold the included masses. The carts also feature threaded inserts for attaching accessories and sensors, and slots for holding a picket fence (such as PASCO part ME-9804) or the included angle indicator.


At each end are two attachment points for connecting a string or spring.

For more information, see the manual packaged with the PAScars.

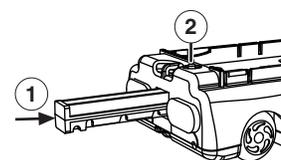

PAScar

- 1: Push plunger in.
- 2: Press button to trigger.

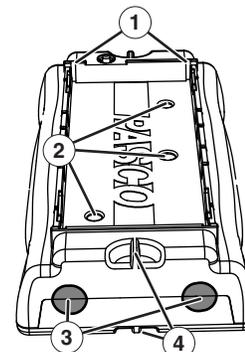

Cart Features

- 1: Slots
- 2: Threaded inserts
- 3: Bumpers
- 4: Attachment points

Metal Carts in ME-5704, ME-5705, and ME-5706


Plunger Cart and Collision Cart These carts run along the track on low-friction wheels.

The plunger cart includes a spring-loaded plunger that launches the cart off of an end stop or another cart. Push the plunger in to the first, second, or third “click” (for a slow, medium, or fast launch); then push up to make the trigger catch. To make the plunger pop out, tap the trigger button with a hard object (such as an included mass bar). When the plunger is not in use, push it all the way in to the third position.

For elastic and inelastic collisions, both carts have hook-and-loop bumpers on one end and magnetic bumpers on the other end. When the hook-and-loop bumpers collide, the carts stick to each other. When the magnetic bumpers collide, the carts repel each other without touching. The magnetic bumpers also repel the magnets in the end stops.


The top of each cart forms a tray designed to hold the included masses. The carts also feature tapped holes for attaching accessories and sensors, and slots for holding a picket fence (such as PASCO part ME-9804) or an angle indicator (PASCO part ME-9495).

For more information, see the manuals packaged with both carts.

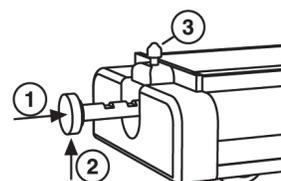

Plunger Cart

- 1: Push plunger in.
- 2: Push up to lock.
- 3: Tap button to trigger.

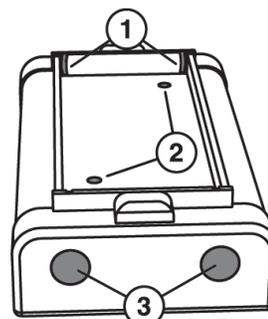


Cart Features

- 1: Slots
- 2: Tapped holes
- 3: Bumpers

Cart Masses


These masses are designed to be placed on the upper tray of a cart.

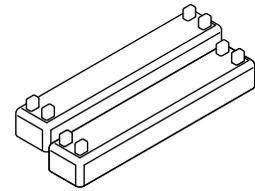

Cart Masses

Pivot Clamp

Attach the pivot clamp to the T-slot of the track (as illustrated) to elevate one end of the track on a vertical rod.

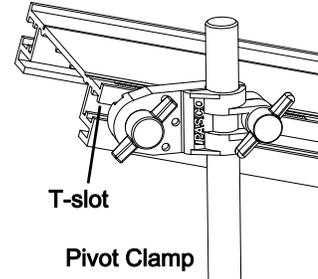

T-slot

Pivot Clamp

Spare Parts

You can order any of the major components of the system using the part numbers listed in the tables for Included Equipment. For an assortment of thumbscrews, square nuts, and other small parts, order the IDS Spares Kit (ME-9823)

For help identifying a part, contact PASCO Technical Support.

Adding Sensors to Your Dynamics System

1. Wireless Force Sensor (one or two)



Wireless Force Acceleration Sensor
PS-3202

We recommend the Wireless Force-Acceleration Sensors because they can ride in the carts without the added friction of a cord. Although a wired force sensor can be used mounted to the end of the track to show impulse, if you want to show the forces between two carts, the wireless force sensors are required.

2. Motion Sensors (2) OR Photogates (2)



Motion Sensor
PS-2103A

Photogate
Heads
ME-9498A
Photogate
Brackets
ME-9806

Picket
Fences
ME-8933

Two Motion Sensors can track the velocity (speed and direction) of both carts throughout a collision for Conservation of Momentum experiments. Photogates have to be placed at just the right position to capture the before and after speeds.

3. An interface for the Motion Sensors or Photogates (550 or 850 Interface)



The 550 is sufficient for most labs, but the 850 will meet every need (particularly for college/university labs).

4. PASCO Capstone Software or SPARKvue

PASCO Capstone Software is best for Windows or Mac in Physics and Engineering laboratories. The site license allows students to use it at home.



Use SPARKvue if you are on a mobile platform: iOS, Android, or Chrome.

Accessories Available Separately

Many additional accessories are available for dynamics systems, such as the Fan Accessory, Friction Block, Magnetic Damping Accessory, and Super Pulley with Clamp. See www.pasco.com for more information.

About the Experiments

These five experiments can be done with the equipment included in the Basic Dynamics Systems and other equipment such as springs, pulleys and mass sets. See each experiment for a specific equipment list.

1. **Conservation of Momentum in Explosions:** Demonstrate conservation of momentum when two carts push off from each other.
2. **Conservation of Momentum in Collisions:** Demonstrate conservation of momentum in elastic and inelastic collisions.
3. **Launch Speed:** Show qualitatively how the final speed of the plunger cart depends on its mass and the initial compression of the plunger spring.
4. **Newton's Second Law :** Show quantitatively how the acceleration of a cart is related its mass and net force.
5. **Acceleration Down an Incline:** Determine the strength of earth's gravity (g) by measuring the acceleration of a cart on an inclined track.

Additional Experiment Suggestions: Many of these suggestions are variations on the above experiments. Some require additional carts.

Conservation of Center of Mass

Set up the track in the configuration shown in Experiment 1 but instead of putting the track directly on the table, place it on a mass bar so that the bar acts as a fulcrum. Position the bar so the carts and track are balanced. First use two carts of equal mass. Trigger the plunger and watch the carts move to the ends of the track. Since the center of mass of the system does not move, the track remains balanced. Repeat this procedure using carts of unequal mass.

Newton's Second Law: Uphill Acceleration

Repeat **Experiment 4** with the track inclined so the pulley is on the high end and the cart accelerates up the incline.

Damped Motion

Incline the track with the end stop at the bottom. Release the cart from a measured distance up the track. Use the magnetic bumpers in the cart and end stop so the cart rebounds. On each rebound, when the cart reaches its peak, record the time and position. Make a graph of amplitude versus time.

Rocket Cart with Balloon

Attach an untied inflated balloon to the cart with the neck of the balloon directed out the back of the cart. Let the air propel the cart.

For the following experiment suggestions, a 2.2 m track (such as model ME-9779) is recommended.

Multiple Inelastic Collisions

Use three carts with the magnets removed from the middle cart. Push the left cart into the middle cart, which in turn will collide with the right cart. The carts will all stick together. Note the resulting final velocity of the carts.

Rocket Staging

Use three or more plunger carts to simulate a rocket expelling fuel. Push the plungers in on each cart and attach the carts together in a line. Use tape to lightly attach the carts to each other. Position the carts at one end of the track. The lead cart represents the rocket and the rest of the carts are fuel. Use a meter stick to release the plungers in succession by striking the plunger-release of each cart, beginning with the last fuel cart (furthest from the rocket cart). As each plunger is released, each cart will separate from the rest, one at a time. Note the final speed of the rocket cart compared to its speed when all the fuel is dumped at once.

Longitudinal Wave

Use six or more carts with magnetic bumpers installed on both ends. Install end stops on the track with the magnetic side toward the center of the track. Start a longitudinal pulse by displacing one of the carts. The carts will rebound off each other and the end stops. Oscillate the end cart to keep a longitudinal wave going down the track.

NOTES

Experiment 1: Conservation of Momentum in Explosions

Required Equipment from Dynamics System

Track with Feet and End Stops	Cart Masses
PASCO Carts (2)	

Other Required Equipment

Suggested Model Number

Mass and Hanger Set	ME-8979
---------------------	---------

Purpose

The purpose of this experiment is to demonstrate conservation of momentum with two carts pushing away from each other.

Theory

When two carts push away from each other (and there is no net force on the system), the total momentum is conserved. If the system is initially at rest, the final momentum of the two carts must be equal in magnitude and opposite in direction to each other so the resulting total momentum of the system is zero:

$$\mathbf{p} = m_1\mathbf{v}_1 - m_2\mathbf{v}_2 = 0$$

Therefore, the ratio of the final speeds of the carts is equal to the ratio of the masses of the carts.

$$\frac{v_1}{v_2} = \frac{m_2}{m_1}$$

To simplify this experiment, the starting point for the carts at rest is chosen so that the two carts will reach the ends of the track simultaneously. The speed, which is the distance divided by the time, can be determined by measuring the distance traveled since the time traveled by each cart is the same.

$$\frac{v_1}{v_2} = \frac{\left(\frac{\Delta x_1}{\Delta t}\right)}{\left(\frac{\Delta x_2}{\Delta t}\right)} = \frac{\Delta x_1}{\Delta x_2}$$

Thus the ratio of the distances is equal to the ratio of the masses:

$$\frac{\Delta x_1}{\Delta x_2} = \frac{m_2}{m_1}$$

Procedure

1. Install the feet on the track and level it. Install one end stop at each end with the magnetic sides of the end stops facing *away* from the carts.

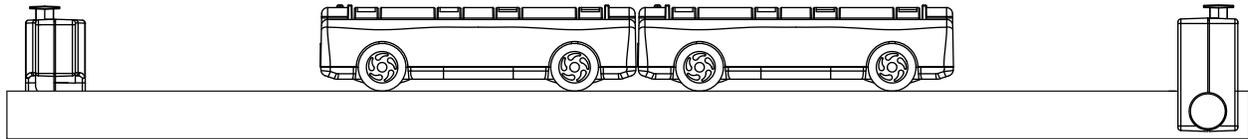


Figure 1.1

2. For each of the cases in Table 1.1, place the two carts against each other with the plunger of one cart latched in the third position and touching the other cart (see Figure 1.1). Push the plunger of the other cart all the way in to its storage position.
3. Tap the plunger release button and watch the two carts move to the ends of the track. Experiment with different starting positions until the two carts reach the ends of the track at the same time. Measure the masses of the carts. Record the masses and the starting position in Table 1.1.

Table 1.1: Results

Additional Mass on Cart 1	Additional Mass on Cart 2	m_1	m_2	Starting Position	x_1	x_2	x_1/x_2	m_2/m_1
0	0							
500 g	0							
1000 g*	0							
500	250							

*Use four 250 g cart masses if possible

Data Analysis

1. For each of the cases, calculate the distances traveled from the starting position to the end of the track. Record the result in Table 1.1.
2. Calculate the ratio of the distances traveled and record in the table.
3. Calculate the ratio of the masses and record in the table.

Questions

1. Does the ratio of the distances equal the ratio of the masses in each of the cases? In other words, is momentum conserved?
2. When carts of unequal masses push away from each other, which cart has more momentum?
3. When the carts of unequal masses push away from each other, which cart has more kinetic energy?
4. Is the starting position dependent on which cart has the plunger? Why?

Experiment 2: Conservation of Momentum in Collisions

Required Equipment from Dynamics System

Track with Feet

PASCO Carts (2)

Other Required Equipment

Paper (for drawing diagrams)

Purpose

The purpose of this experiment is to qualitatively explore conservation of momentum for elastic and inelastic collisions.

Theory

When two carts collide with each other, the total momentum of both carts is conserved regardless of the type of collision.

An elastic collision is one in which the carts bounce off each other with no loss of kinetic energy. In this experiment, magnetic bumpers are used to minimize the energy losses due to friction during the collision. In reality, this “elastic” collision is slightly inelastic.

A completely inelastic collision is one in which the carts hit and stick to each other. In this experiment, this is accomplished with the hook-and-loop bumpers on the carts.

Part I: Elastic Collisions

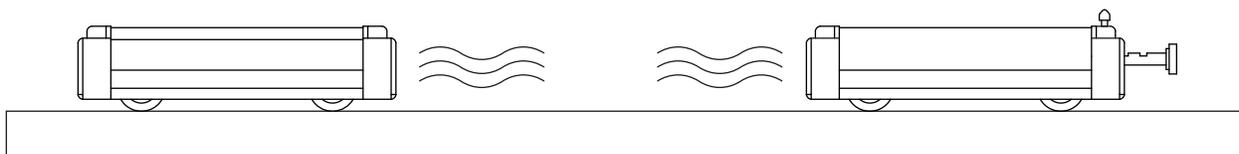


Figure 2.1

1. Install the feet on the track and level it.
2. Orient the two carts on the track so their magnetic bumpers are toward each other.
3. Test cases A1 through A3 and B1 through B3 described below. Draw two diagrams (one for before the collision and one for after the collision) for each case. In every diagram, show a velocity vector for each cart with a length that approximately represents the relative speed of the cart.

A. Carts with Equal Mass

Case A1: Place one cart at rest in the middle of the track. Give the other cart an initial velocity toward the cart at rest.

Case A2: Start the carts with one at each end of the track. Give each cart approximately the same velocity toward each other.

Case A3: Start both carts at one end of the track. Give the first cart a slow velocity and the second cart a faster velocity so that the second cart catches the first cart.

B. Carts with Unequal Mass

Put two mass bars in one of the carts so that the mass of one cart is approximately three times the mass (3M) of the other cart (1M).

Case B1: Place the 3M cart at rest in the middle of the track. Give the other cart an initial velocity toward the cart at rest.

Case B2: Place the 1M cart at rest in the middle of the track. Give the 3M cart an initial velocity toward the cart at rest.

Case B3: Start the carts with one at each end of the track. Give each cart approximately the same velocity toward each other.

Case B4: Start both carts at one end of the track. Give the first cart a slow velocity and the second cart a faster velocity so that the second cart catches the first cart. Do this for both cases: with the 1M cart first and then for the 3M cart first.

Part II: Completely Inelastic Collisions

1. Orient the two carts so their hook-and-loop bumpers are toward each other. Push the plunger in all the way so it will not interfere with the collision.
2. Repeat test cases A1 through A3 and B1 through B3 and draw diagrams for each case.

Questions

1. When two carts having the same mass and the same speed collide and stick together, they stop. Is momentum conserved?
2. When two carts having the same mass and the same speed collide and bounce off of each other elastically, what is the final total momentum of the carts?

Experiment 3: Launch Speed

Required Equipment from Dynamics System

Track with Feet and End Stops	Cart Masses
PASCO Cart	Pivot Clamp

Other Required Equipment

Suggested Model Number

Mass and Hanger Set	ME-8979
---------------------	---------

Purpose

In this experiment, you will qualitatively demonstrate how the final speed of the plunger cart depends on its mass and the initial compression of the plunger spring.

Procedure

1. Install the feet on the track and level it.
2. Install an end stop at each end of the track.
3. Push the plunger of the cart into the first, second, or third position. Place the cart on the track with the plunger against one end stop. Tap the trigger button to launch the cart.
4. **Varying Spring Compression:** Repeat step 3 three times. Perform the first trial with the spring plunger cocked to the first possible position (the least compression) and then do two more trials increasing the force applied to the cart by increasing the compression of the spring plunger.
5. **Varying Mass:** For these trials, always cock the spring plunger to the maximum. Observe the relative accelerations of the cart alone and with 250 g, 500g, 750 g, and 1000 g of added mass.

Questions

1. When does the spring apply force to the cart? When is the net force on the cart (ignoring friction) zero?
2. Does the final speed of the cart increase or decrease with increasing initial spring compression?
3. Does the final speed of the cart increase or decrease with increasing cart mass?

NOTES

Experiment 4: Newton's Second Law

Required Equipment from Dynamics System

Track with Feet and End Stop	Cart Masses
PASCO Cart	

Other Required Equipment

Other Required Equipment	Suggested Model Number
Super Pulley with Clamp	ME-9448
Stopwatch	ME-1234
Mass and Hanger Set	ME-8979
Mass balance	SE-8723
String (about 2 m)	SE-8050

Purpose

In this experiment, you will verify Newton's Second Law, $F = ma$.

Theory

According to Newton's Second Law, $F = ma$, where F is the net force acting on the object of mass m , and a is the resulting acceleration of the object.

For a cart of mass m_1 on a horizontal track with a string attached over a pulley to a hanging mass m_2 (see Figure 4.1), the net force F on the entire system (cart and hanging mass) is the weight of hanging mass, $F = m_2g$, (assuming that friction is negligible).

According to Newton's Second Law, this net force should be equal to ma , where m is the total mass that is being accelerated, which in this case is $m_1 + m_2$. You will check to see if $m_2g = (m_1 + m_2)a$ as predicted by theory.

To determine the acceleration, you will release the cart from rest and measure the time (t) for it to travel a certain distance (d). Since $d = (1/2)at^2$, the acceleration can be calculated using $a = 2d/t^2$.

Procedure

1. Install the feet on the track and level it.
2. Install one end stop on the track near one end with the magnets facing away from the track.
3. Measure the mass of the cart and record it in Table 4.1.

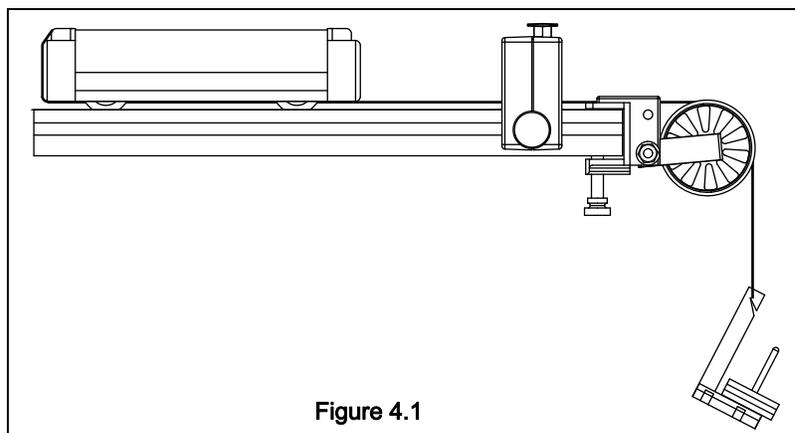


Figure 4.1

4. Attach the pulley and end stop to the track as shown in Figure 4.1. Place the cart on the track. Tie a string to the lower tie point of the cart. Tie a mass hanger on the other end of the string. Run the string under the end stop and over the pulley. Adjust the pulley so that the string runs parallel to the track. The string must be just long enough so the cart reaches the end stop before the mass hanger reaches the floor.

- Pull the cart back until the mass hanger reaches the pulley. Record this initial release position in Table 4.1. This will be the release position for all the trials. Make a test run to determine how much mass is required on the mass hanger so that the cart takes about 2 seconds to complete the run. Because of reaction time, too short of a total time will cause too much error. However, if the cart moves too slowly, friction causes too much error. Record the hanging mass in Table 4.1.
- Place the cart against the end stop on the pulley end of the track and record the final position of the cart in Table 4.1.
- Pull the cart back to the initial release position. Release it and time how long it takes to reach the end stop. Record the time in Table 4.1.
- Measure the time at least 5 times with the same mass and record these values in Table 4.1.

Table 4.1: Experimental Period

Initial release position = _____							
Final position = _____							
Distance traveled (d) = _____							
Cart Mass	Hanging Mass	Time					Average Time
		Trial 1	Trial 2	Trial 3	Trial 4	Trial 5	

- Add a mass bar to the cart and repeat the procedure.

Data Analysis

- Calculate the average times and record them in Table 4.1.
- Record the distance traveled (from initial to final position) in Table 4.1.
- Calculate the accelerations and record them Table 4.2.
- For each case, calculate $(m_1 + m_2)a$ and record in Table 4.2.
- For each case, calculate the net force, F_{NET} , and record in Table 4.2.
- For each case, calculate the percent difference between F_{NET} and $(m_1 + m_2)a$ and record in Table 4.2.

Table 4.2

Cart Mass	Acceleration	$(m_1 + m_2)a$	$F_{\text{NET}} = m_2g$	% Difference

Questions

- Did the results of this experiment verify that $F = ma$?
- Why must the mass in $F = ma$ include the hanging mass as well as the mass of the cart?

Experiment 5: Acceleration Down an Incline

Required Equipment from Dynamics System

Track with End Stop	Pivot Clamp
PASCO Cart	

Other Required Equipment

Other Required Equipment	Suggested Model Number
Base and support rod	ME-9355
Stopwatch	ME-1234
Graph paper	

Purpose

In this experiment, you will investigate how the acceleration of a cart rolling down an inclined track depends on the angle of incline. From your data, you will calculate the acceleration of an object in free-fall.

Theory

A cart of mass m on an incline will roll down the incline as it is pulled by gravity. The force of gravity (mg) is straight down as shown in Figure 5.1. The component of that is parallel to the inclined surface is $mg \sin \theta$.

To determine the acceleration, you will release the cart from rest and measure the time (t) for it to travel a certain distance (d). Since $d = (1/2)at^2$, the acceleration can be calculated using $a = 2d/t^2$.

A plot of a versus $\sin \theta$ will be a straight line with a slope equal to the acceleration of an object in free-fall, g .

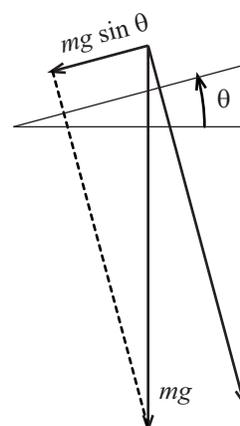


Figure 5.1

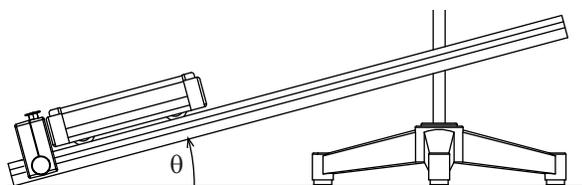


Figure 5.2

Procedure

1. Set up the track as shown in Figure 5.2 with a pivot clamp and support stand. Elevate the end of the track by about 10 cm.
2. Set the cart on the track against the end stop and record this final position in Table 5.1. (Use the non-magnetic end of the cart so it touches the end stop.)
3. Pull the cart up to the top of the track and record the initial position where the cart will be released from rest.
4. Release the cart from rest and use the stopwatch to time how long it takes the cart to reach the end stop. The person who releases the cart should

also operate the stopwatch. Repeat this measurement 10 times (with different people doing the timing). Record all the values in Table 4.1.

- Lower the end of the track by 1 cm and repeat step 4. Use the same release position.
- Repeat step 4 for a total of 7 angles, lowering the end of the track by 1 cm for each new angle.

Table 5.1

		Initial release position = _____ Final position = _____ Distance traveled (d) = _____						
		Height of Track						
		10 cm	9 cm	8 cm	7 cm	6 cm	5 cm	4 cm
Time	Trial 1							
	Trial 2							
	Trial 3							
	Trial 4							
	Trial 5							
	Trial 6							
	Trial 7							
	Trial 8							
	Trial 9							
	Trial 10							
	Average							

Data Analysis

- Calculate the average time for each angle and record it in Table 5.1.
- Calculate the distance traveled, d , from the initial to the final position.
- Use the distance traveled and average time to calculate the acceleration for each angle and record it in in Table 5.2.

Table 5.2

Height	Acceleration	$\sin \theta$
10 cm		
9 cm		
8 cm		
7 cm		
6 cm		
5 cm		
4 cm		

4. Measure the hypotenuse of the triangle formed by the track and use this to calculate $\sin \theta$ for each angle.
 - Hypotenuse = _____
5. Plot acceleration versus $\sin \theta$. Draw the best-fit straight line and calculate its slope. Calculate the percent difference between the slope and $g = 9.8 \text{ m/s}^2$.
 - slope = _____
 - % difference = _____

Questions

1. Does your reaction time in operating the stopwatch cause a greater percentage error at higher or lower track angles?
2. How will doubling the mass of the cart affect the results? Try it.

NOTES

Technical Support

For assistance with any PASCO product, contact PASCO at:

Address: PASCO scientific
10101 Foothills Blvd.
Roseville, CA 95747-7100

Phone: 916-462-8384 (worldwide)
800-772-8700 (U.S.)

Web: www.pasco.com

Email: support@pasco.com

Limited Warranty

For a description of the product warranty, see the PASCO catalog.

Copyright

This PASCO scientific *Instruction Manual* is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

Trademarks

PASCO and PASCO scientific are trademarks or registered trademarks of PASCO scientific, in the United States and/or in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of, their respective owners. For more information visit www.pasco.com/legal.

