



**Instruction Sheet
for the PASCO
Model CI-6692**

IDS MOUNT ACCESSORY

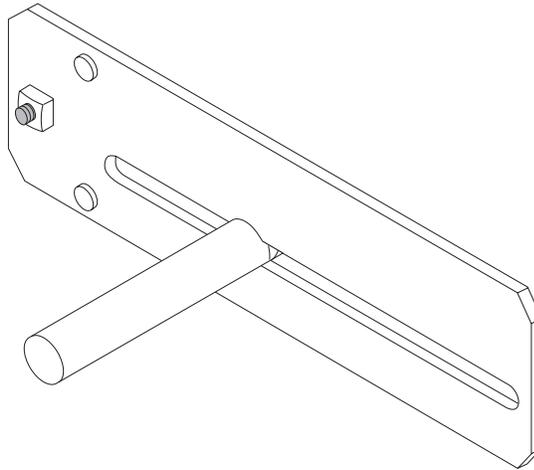


The PASCO CI-6692 IDS Mount Accessory is designed for use with the PASCO CI-6538 Rotary Motion Sensor or the PASCO CI-6625 Rotary Motion Sensor for ULI. This bracket allows the Rotary Motion Sensor to be easily attached to the PASCO Introductory Dynamics System tracks (1.2 meter ME-9435A or 2.2 meter ME-9458).

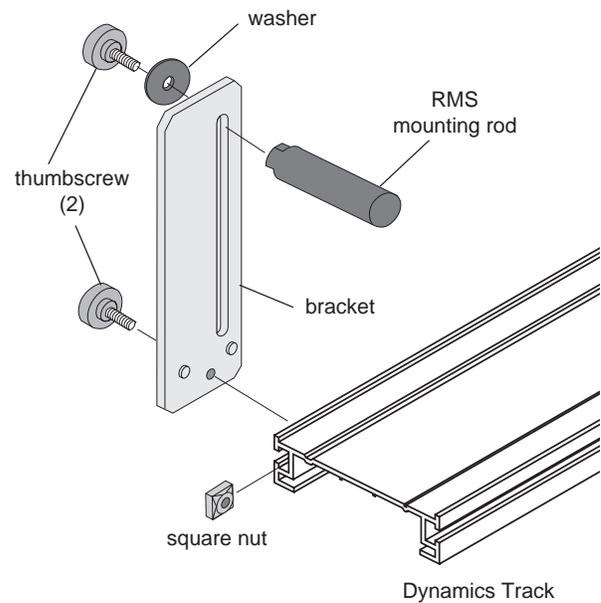

Attaching the Bracket to the Track

© 1996 PASCO scientific

This instruction sheet written by: Jon Hanks

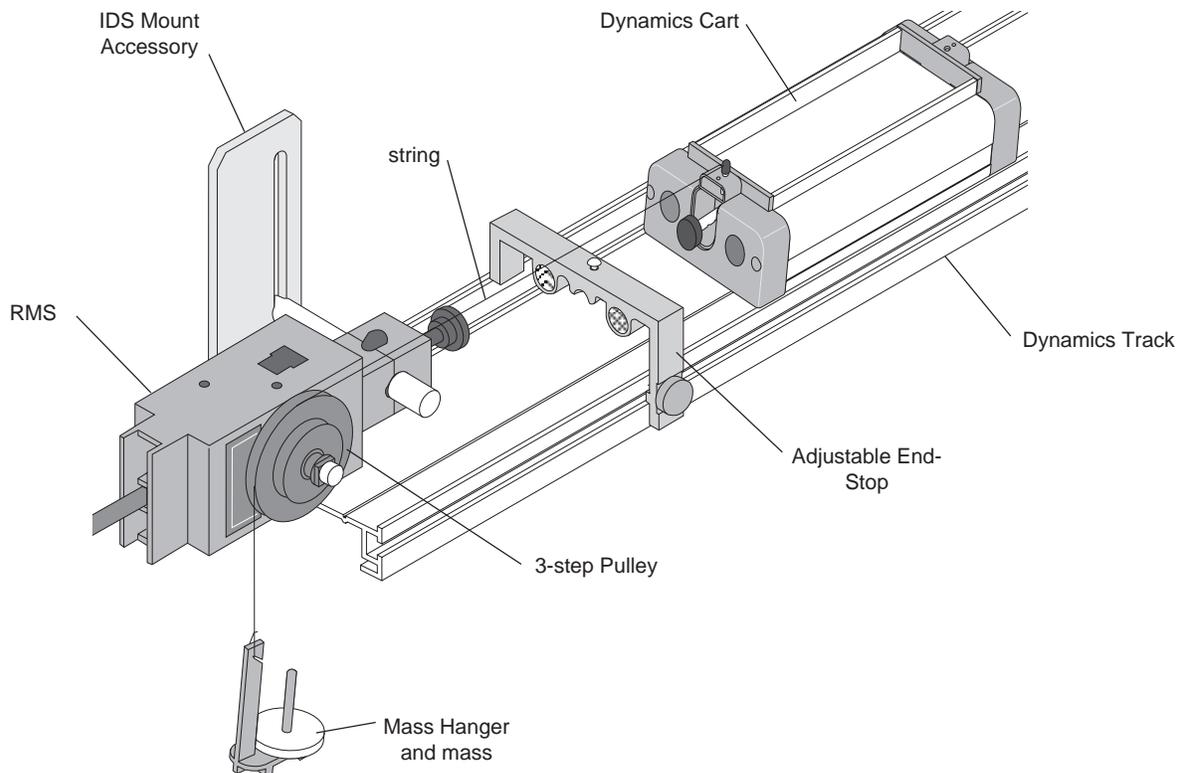

Equipment Setup Using RMS Like a Smart Pulley

Limited Warranty

PASCO scientific warrants this product to be free from defects in materials and workmanship for a period of one year from the date of shipment to the customer. PASCO will repair or replace, at its option, any part of the product which is deemed to be defective in material or workmanship. This warranty does not cover damage to the product caused by abuse or improper use. Determination of whether a product failure is the result of a manufacturing defect or improper use by the customer shall be made solely by PASCO scientific. Responsibility for the return of equipment for warranty repair belongs to the customer. Equipment must be properly packed to prevent damage and shipped postage or freight prepaid. (Damage caused by improper packing of the equipment for return shipment will not be covered by the warranty.) Shipping costs for returning the equipment, after repair, will be paid by PASCO scientific.

To Reach PASCO

For Technical Support call us at 1-800-772-8700 (toll-free within the U.S.) or (916) 786-3800.

email: techsupp@PASCO.com

Tech support fax: (916) 786-3292

Web: <http://www.pasco.com>