


AirLink Replacement Battery

PS-2566

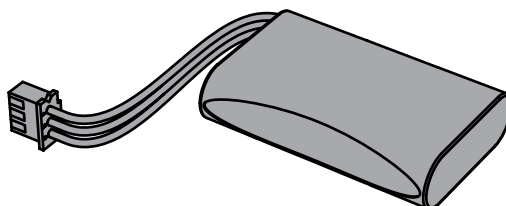

Introduction

PASCO model PS-2566 is a replacement battery for the PASPORT AirLink SI.

To replace the battery, you will need to open your AirLink and connect the new battery to the circuit board. If you would prefer to send your AirLink back to PASCO to have the battery replaced, contact PASCO by phone or email (see Technical Support on page 3).

What You Will Need

To replace the battery, you will need the following items.


- AirLink Replacement Battery (PASCO model PS-2566)
- AirLink SI (PASCO model PS-2005)
- T5 Torx driver

Replacing the Battery


Important: Before proceeding, take precautions to avoid static discharge, which may damage the AirLink. Touch something metal and grounded before handling the AirLink circuit board.

Opening the AirLink

1. Using a T5 Torx driver, remove the two screws from the bottom half of the AirLink enclosure.

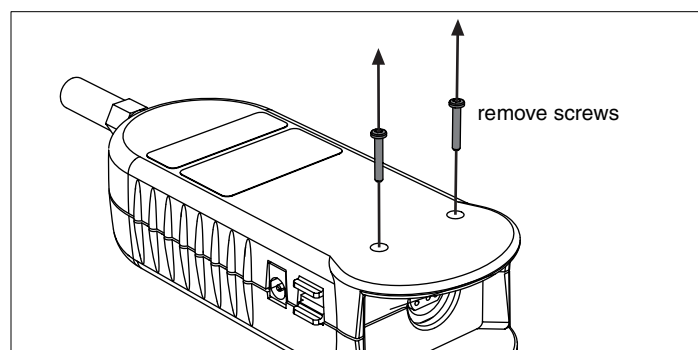


2. Firmly pull apart the halves of the enclosure. As you open the AirLink, the circuit board will separate from the top half of the enclosure.

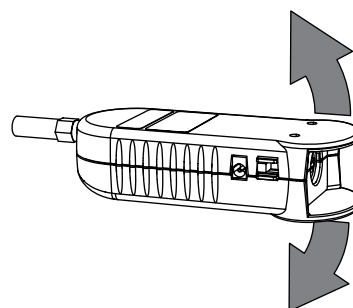


Removing the Old Battery


3. Disconnect the battery plug by pulling it directly away from the circuit board.

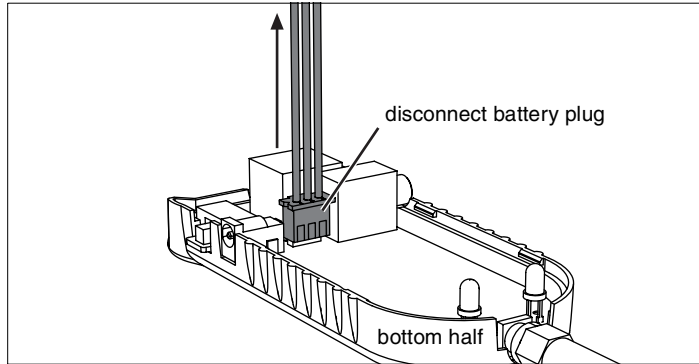

4. Replace the circuit board in the bottom half of the enclosure and set both parts aside.
5. Using a T5 Torx driver, remove the two screws securing the battery cover.

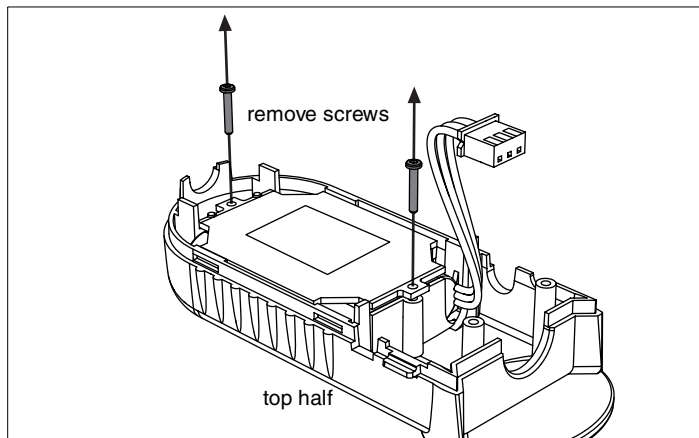


6. Remove the battery cover.
7. Remove the old battery.

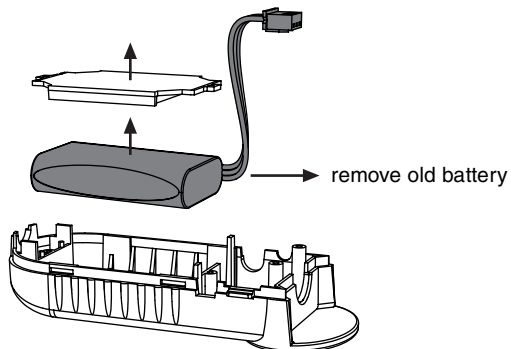


Installing the New Battery

8. Place the new battery in the battery compartment as illustrated.

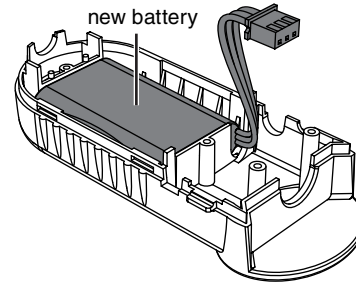

Important: Do not over tighten the screws. Tighten them only enough to hold the battery cover in place.


9. Replace the battery cover. Be sure to align the notches with the posts as illustrated.

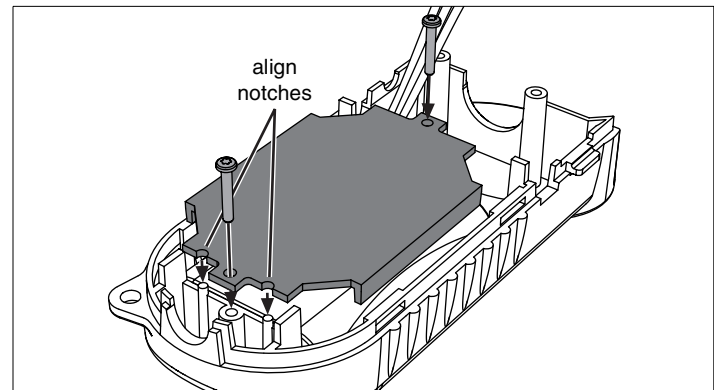

10. Secure the battery cover with two screws.
11. Observe the correct orientation of the battery plug and wires illustrated below.


Important: Do not connect the battery backwards. The battery may be damaged if it is connected incorrectly.

12. Connect the battery plug to the circuit board as illustrated.

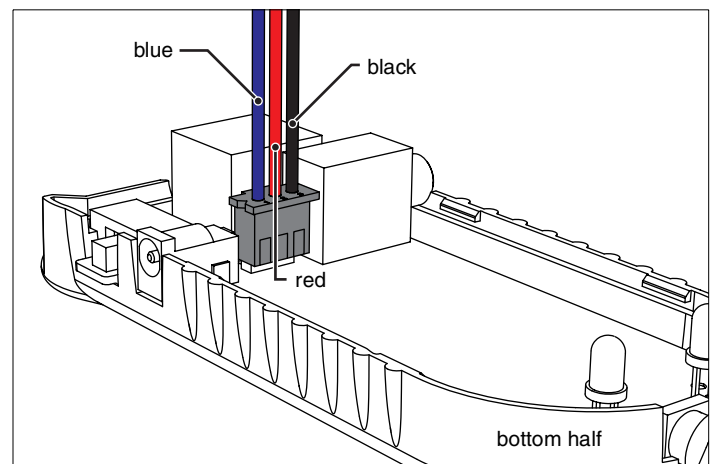

Reassembling the AirLink


13. Place the bottom half of the AirLink (containing the circuit board) flat on a tabletop.

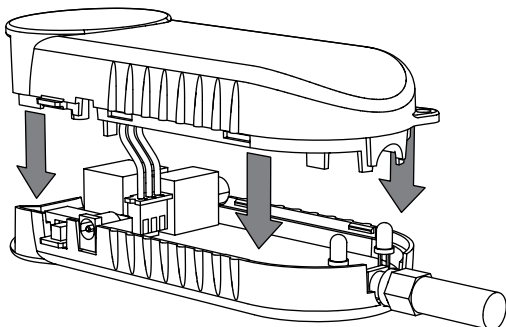

14. Align the halves of the AirLink and press firmly to snap them together.

Important: Do not over tighten the screws. Tighten them only enough to seat them in place.

15. Replace the two screws in the enclosure.

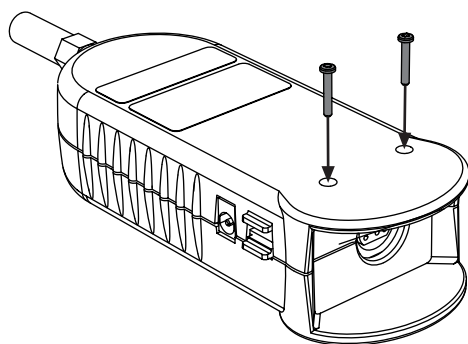

Charging the New Battery

16. Set the switch on the AirLink to Off.
17. Connect the AC adapter (included with the AirLink) to the DC Power input of the AirLink.
18. Connect the AC adapter to a wall socket.

A green LED in the AirLink blinks as the battery charges. When the battery is fully charged (after about four hours), the LED glows steadily.

Technical Support

If you have questions or if you need replacement parts, please contact PASCO technical support.

Address: PASCO scientific
10101 Foothills Blvd.
Roseville, CA 95747-7100

Phone: 916-786-3800 (worldwide)
800-772-8700 (U.S.)

Fax: (916) 786-7565

Web: www.pasco.com

Email: support@pasco.com

Limited Warranty For a description of the product warranty, see the PASCO catalog.

Copyright The PASCO scientific 012-10290A *AirLink Replacement Battery Instruction Sheet* is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

Trademarks PASCO, PASCO scientific, and PASPORT are trademarks or registered trademarks of PASCO scientific, in the United States and/or in other countries. Torx is a registered trademark of Acument Intellectual Properties, LLC. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of, their respective owners. For more information visit www.pasco.com/legal.